

Ebola Pseudoviral Particles – Zaire Ebolavirus (Makona 2014 Strain)

CATALOG NUMBER: PsLv-EBOV-M2014

Description

Ebola virus enters host cells via its surface glycoprotein (GP), which binds to specific receptors on the cell membrane and facilitates membrane fusion.

eEnzyme's PsLv-EBOV-M2014 pseudoviral particles are replication-deficient lentiviruses pseudotyped with the GP from Sudan ebolavirus, strain H.sapiens-wt/GIN2014/Makona-Coyah-955 (GenBank Accession No. AKG65250). These particles encode firefly luciferase (FLuc) as a reporter, enabling quantification of viral entry through luminescence. This system isolates the entry step from the rest of the viral life cycle, making it ideal for targeted assays.

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


Figure 1. Illustration of the replication-deficient lentivirus pseudotyped with Ebola glycoproteins

Applications

Optimized for use with eEnzyme's Firefly Luciferase Assay Kit (Catalog # CA-L165), these pseudoviruses are ideal for:

- Screening inhibitors of Ebola virus entry and GP-mediated fusion
- Evaluating neutralizing antibodies against Zaire ebolavirus
- High-throughput infection assays in 96-, 384-, and 1536-well formats

Key Features

- **Robust Signal:** High signal-to-noise ratio for sensitive detection
- **HTS-Compatible:** Validated for multi-well formats
- **Convenient Format:** Ready-to-use for luciferase-based entry assays

Product Contents

- Volume: 1 mL or 5 mL per vial
- Format: Lentivirus-based pseudoviral particles
- Titer: $>1.0 \times 10^7$ particles/mL

Storage & Stability:

- Store immediately at -70°C upon receipt



- Thaw rapidly at room temperature (<30 min, no shaking); use immediately
- Do not aliquot or refreeze
- Shelf life: 6 months from shipping date when stored at -70°C

Assay Protocol

Cell Infection

1. Seed HEK293 cells:
 - 96-well: 20K cells/well in 50 μ L DMEM + 10% HyClone™ FetalClone™ II Serum (no antibiotics)
 - 384-well: 5K cells/well in 15 μ L
2. Incubate overnight for adherence
3. Replace media and add pseudoviral particles:
 - 96-well: 50 μ L/well
 - 384-well: 12.5 μ L/well
 - Centrifuge at 700 rpm for 15 min at 4°C
4. Incubate for 2 hrs at 37°C
5. Add fresh DMEM + 10% serum (same volume as above)
6. Incubate for 48-72 hrs at 37°C

Luciferase Readout

1. Without removing supernatant, add luciferase assay working solution:
 - 96-well: 100 μ L/well
 - 384-well: 25 μ L/well
 - Use eEnzyme's Firefly Luciferase Assay Kit (Catalog # CA-L165)
2. Measure luminescence using a microplate reader (e.g., BioTek Synergy 2, gain 255)

Representative Data

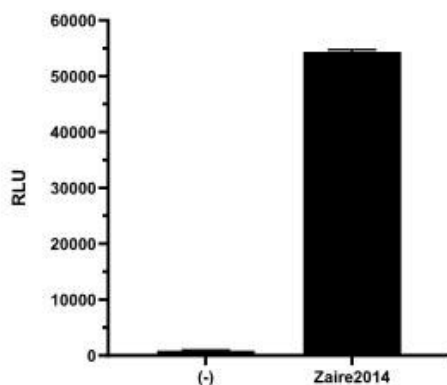


Figure 2: Assay: HEK293 cells infected with PsLv-EBOV-M2014 pseudovirus

Readout: Relative Luminescence Units (RLU)

(-): Negative control using lentivirus without envelope protein

Zaire2014: Ebola GP pseudovirus (PsLv-EBOV-M2014)

